



STEM CELL LABORATORY (STCL)



DOCUMENT NUMBER: STCL-PROC-018

DOCUMENT TITLE:

Correct Connection of Pre-system Filter and Tubing Set-CliniMACS

DOCUMENT NOTES:

FACT # 6D.700.

Document Information

Revision: 03

Vault: STCL-Processing-rel

Status: Release

Document Type: STCL-Processing

Date Information

Creation Date: 03 May 2011

Release Date: 21 Jan 2013

Effective Date: 21 Jan 2013

Expiration Date:

Control Information

Author: WATE02

Owner: WATE02

Previous Number: 6D.700 \JA1\


Change Number: FORM-CCR-119

STCL-PROC-018 CORRECT CONNECTION OF PRE-SYSTEM FILTER AND TUBING SET


1. Correct spiking depth into the pre-system filter

When inserting the tubing set spike into the pre-system filter port, please ensure that the aperture of the spike is completely covered by the port. The connection is safe when approx. 5-8 mm of the spike below the aperture is inserted (white “ring” caused by direct contact of spike and port must be visible)

Spike too deep inserted

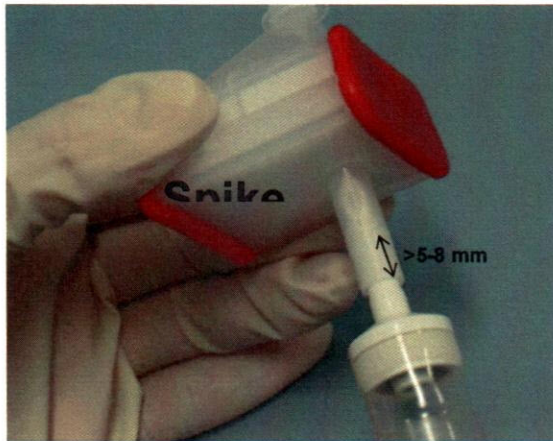


Correct spike insertion

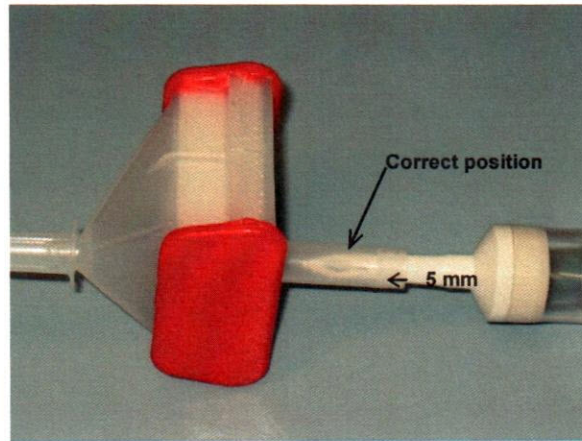


2. How to hold the tubing set spike and filter during insertion

For a correct connection, insert the spike in a half-turn by holding the drip chamber in one hand and the pre-system filter in the other hand. It is important not to hold the tubing set spike behind the drip chamber and the pre-system filter at the rear spike as both hands have to be close together during insertion to avoid breakages.

**Hands too far away from each other
will increase breakage risk**

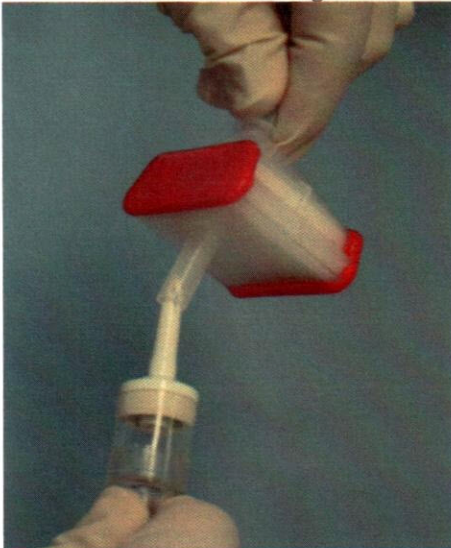

**Hands close together ensure correct insertion
without risk of breakage**

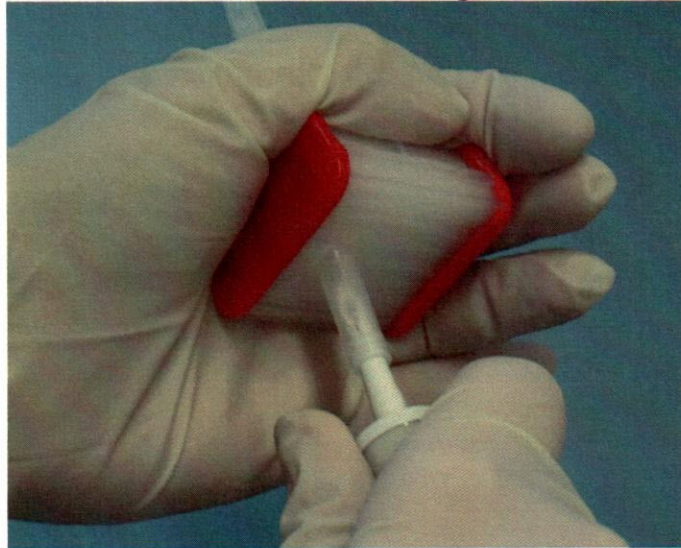

Signature Manifest**Document Number:** STCL-PROC-018**Revision:** 03**Title:** Correct Connection of Pre-system Filter and Tubing Set-CliniMACS**STCL-PROC-018 Pre-System Filter****Author Approval**

Name/Signature	Title	Date	Meaning/Reason
Barbara Waters-Pick (WATE02)		18 Dec 2012, 09:19:30 PM	Approved

Manager Approval

Name/Signature	Title	Date	Meaning/Reason
Barbara Waters-Pick (WATE02)		18 Dec 2012, 09:19:53 PM	Approved

Medical Director Approval

Name/Signature	Title	Date	Meaning/Reason
Joanne Kurtzberg (KURTZ001)		18 Dec 2012, 11:06:08 PM	Approved

QA Approval

Name/Signature	Title	Date	Meaning/Reason
Linda Sledge (SLEDG006)		19 Dec 2012, 08:13:30 AM	Approved

Document Release

Name/Signature	Title	Date	Meaning/Reason
Sandy Mulligan (MULLI026)		18 Jan 2013, 10:04:05 AM	Approved

Notification

Name/Signature	Title	Date	Meaning/Reason
Ann Kaestner (KAEST002)		18 Jan 2013, 10:04:05 AM	Email Sent
Barbara Waters-Pick (WATE02)		18 Jan 2013, 10:04:05 AM	Email Sent